

First time users: Configuring Coherent Connect for start up (UV laser)

- The 4 OBIS model lasers are set to turn on along with the A3 power button. You do not need to start these up, or view them in CC
- Here are the steps to configure **only** the UV (Genesis G4) laser to open up in CC.

**Note: for troubleshooting missing laser signals, refer to Help Files document. This will instruct you to use a different log in that is configured for the 5 CC laser windows to show up.*

At your first visit:

1. Start with A3 powered off: You will start it up after configuring CC, and check by bead QC that lasers turned on OK
2. Log in to PC (AD log in), open Coherent Connect.
3. Go to Laser Options (a) menu
4. Change selections (b) and (c) according to the image on the right.
5. Click OK (d), exit CC
6. See [A3 Quick Guide](#) (printed copy on desk top): Follow along and Start up the A3 according to the guide.
7. After UV laser is started up in CC, perform bead QC as per the A3 Quick Guide to verify CC configuration is working correctly (i.e. all lasers powered on, no missing signals for CST beads).

