



Troubleshooting Red Laser Off: Logging into BdAdmin Account

1. If you are already logged into the computer, start by logging out, and logging into the BdAdmin account under local domain.

2. To sign into local domain BdAdmin Account type in FCCF-SYMPHONY\BdAdmin for the username and BDIS#2\$\$ for the password.

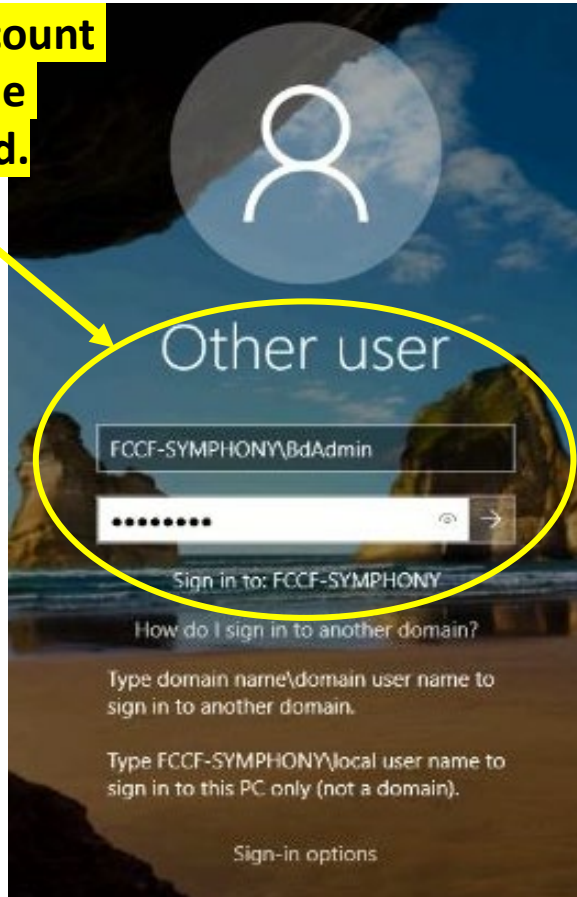

3. Once you are logged on, open Coherent Connect, then wait until you see all 5 laser windows. Turn on the UV and Red laser in the two-step process (1) turning the slide bar up (2) pressing laser START. See that the red dial moves, showing laser is ON.

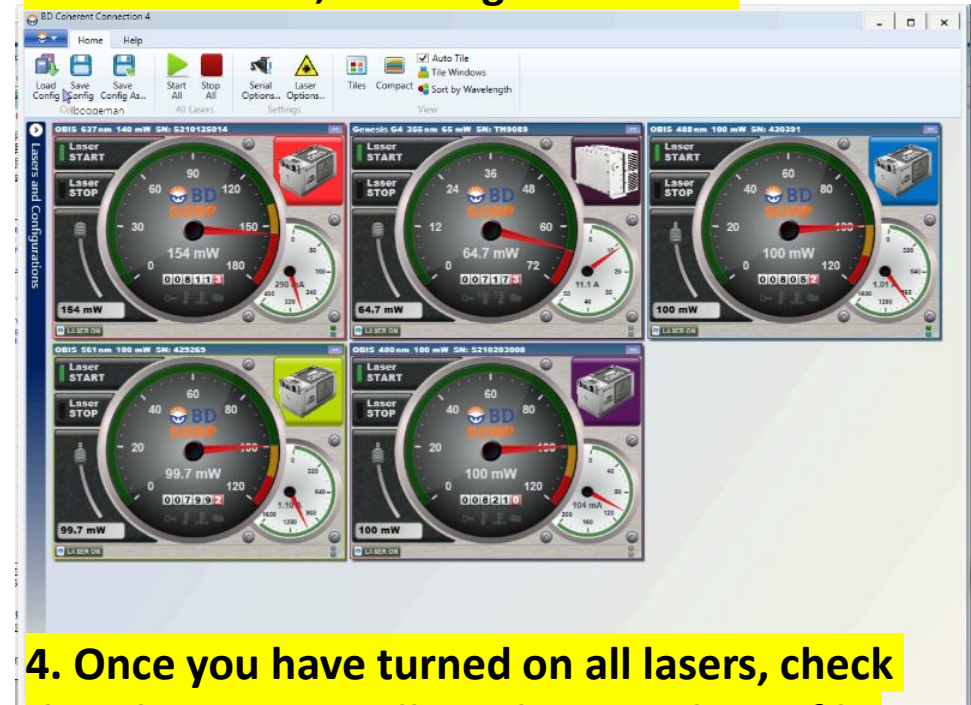

4. Once you have turned on all lasers, check that they are visually on by recording a file of the CST beads histograms in Diva.